

5.90" x 5.90", 1.18" Focal Length, Fresnel Lens



Aspherically Contoured Fresnel Lenses

Stock #13-459 **5 In Stock**

⊖ 1 ⊕ £95⁰⁰

ADD TO CART

| Volume Pricing | |
|----------------|-------------------------------|
| Qty 1-10 | £95.20 each |
| Qty 11-49 | £85.60 each |
| Need More? | Request Quote |

ⓘ Prices shown are exclusive of VAT/local taxes

Product Downloads

General

Fresnel Lens **Type:**

Physical & Mechanical Properties

0.08 **Center Thickness CT (inches):**

±0.05 **Dimensional Tolerance (inches):**

5.9 x 5.9 **Dimensions (inches):**

150.0 x 150.0 **Dimensions (mm):**

±40 **Thickness Tolerance (%):**

Optical Properties

30.00 **Effective Focal Length EFL (mm):**

Acrylic **Substrate:**

Uncoated **Coating:**

400 - 1100 **Wavelength Range (nm):**

1.18 **Effective Focal Length EFL (inches):**

125.00 **Groove Density (grooves/inch):**

1.49 **Index of Refraction (n_d):**

85 (Typical) **Transmission (%):**

Environmental & Durability Factors

80 (Maximum) **Operating Temperature (°C):**

Regulatory Compliance

[Compliant](#) **RoHS 2015:**

[Compliant](#) **Reach 219:**

[View](#) **Certificate of Conformance:**

Product Details

- Positive Focal Length
- Square and Rectangular Options
- Ideal for Light Gathering Applications

Fresnel Lenses replace the curved surface of a conventional lens with a series of concentric grooves, molded into the surface of a thin, lightweight plastic sheet. The grooves act as individual refracting surfaces, like tiny prisms when viewed in cross section, bending parallel rays in a very close approximation to a common focal length. Because the lens is thin, very little light is lost by absorption. Fresnel Lenses are a compromise between efficiency and image quality. High groove density allows higher quality images, while low groove density yields better efficiency (as needed in light gathering applications). In infinite conjugate systems, the grooved side of the lens should face the longer conjugate.

Fresnel lenses are most often used in light gathering applications, such as condenser systems or emitter/detector setups. Fresnel lenses can also be used as magnifiers or projection lenses; however, due to the high level of distortion, this is not recommended.

Technical Information

