

[See all 62 Products in Family](#)

600 Grooves, 12.7mm Square, 1250nm Ruled Diffraction Grating



Reflective Ruled Diffraction Gratings

Stock #43-746 **13 In Stock**

⊖ 1 ⊕ £76.⁰⁰

ADD TO CART

Volume Pricing

Qty 1-9	£76.00 each
Qty 10-24	£68.40 each
Need More?	Request Quote

ⓘ Prices shown are exclusive of VAT/local taxes

Product Downloads

General

Reflective Diffraction Grating **Type:**

Physical & Mechanical Properties

12.7 x 12.7 ±0.5 **Dimensions (mm):**

90 **Clear Aperture (%):**

Ruled Grating	Construction:
Parallel to Short Dimension	Direction of Grooves:
12.70	Length (mm):
6.00 ±0.5	Thickness (mm):
12.70	Width (mm):
±0.5	Alignment of Grooves to Edge (°):

Optical Properties	
600	Groove Density (grooves/mm):
1250	Blaze Wavelength (nm):
22.02	Blaze Angle (°):
Bare Aluminum	Coating:
Float Glass	Substrate: <input type="checkbox"/>
88	Absolute Peak Efficiency, Typical (%):
>83	Peak Efficiency Avg, Typical (%):

Regulatory Compliance	
Compliant	RoHS 2015:
View	Certificate of Conformance:
Compliant	Reach 247:

Product Details

- Up to 80% Efficiency
- Bare Aluminum Coating
- Blaze Wavelength Options from 250 to 1600nm

Reflective Ruled Diffraction Gratings are optical components that separate (diffract) white (polychromatic) light into its component wavelengths. Theoretical resolving power is directly proportional to the number of grooves per millimeter ruled into the master grating. The blaze angle and groove spacing determine the specific wavelength and spectral region where the grating has its maximum efficiency. Reflective Ruled Diffraction Gratings are high quality, instrument-grade gratings that are ideal when high efficiency is the primary concern in devices such as monochromators and spectrometers. Each grating is fabricated from a highly accurate master grating that is copied many times. The duplication process is described below for replicated gratings.

Diffraction ruled gratings are used in a variety of monochromators for research, student and industry use. Almost all commercially available spectrophotometers (ultraviolet, visible, infrared, fluorescence, Raman, atomic absorption) utilize diffraction gratings to select specific wavelengths or scan over a wavelength interval. Generally, replicated ruled gratings should be used when high peak efficiency and throughput is required. Replicated holographic gratings should be used when minimum stray light is critical and high resolution is needed. Note: Damage thresholds for both ruled and holographic gratings are 350 milli-joules/cm² for pulsed lasers and 40 Watts/cm² for CW lasers.


Handling Gratings: Gratings require special handling, making them prone to fingerprints and aerosols. Gratings should only be handled by the edges. Before attempting to clean a grating, please [contact us](#).

Manufacturing of Ruled Gratings

Ruled diffraction gratings are produced by ruling a series of closely spaced, straight parallel grooves into an optically flat aluminum coated substrate, known as the "master grating." Precise, interferometrically-controlled ruling engines utilize a very fine diamond tool to form a sawtooth-shaped groove profile at a given angle (commonly referred to as the blaze angle) on the surface of a prepared substrate. The replication process begins with the surface contour of a ruled master grating being vacuum deposition-coated with an extremely thin separation layer. An aluminum coating is then deposited on top of this separation layer. Then, an epoxy-coated flat glass substrate is placed on top of the layer-covered master, duplicating the grooved surface. The combination is cured and the process is finished when the replicated grating is separated from the master grating.

Technical Information

Typical Efficiency Curves for Ruled Gratings
Optimized (Blaze) Wavelengths from 1000-1600 nm

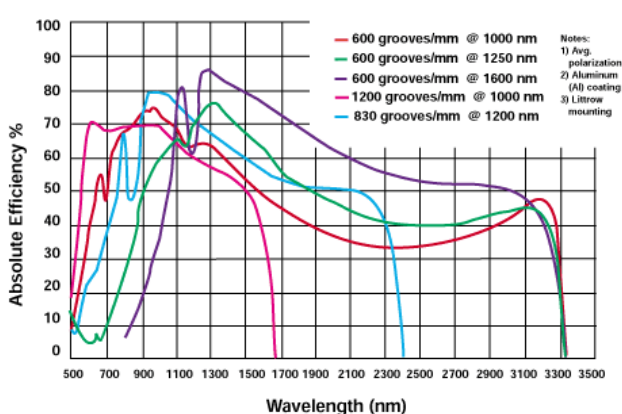

Special Handling

These optics require special handling to avoid damage and ensure long-term performance. Proper handling, cleaning, and storage are essential to maintain optical quality. Explore our [Optics Cleaning Resources](#) for step-by-step guides and best practices. For personalized assistance, [Email us](#) or [Chat](#) with our technical support team.



Component Handling Tools

;